

## Uniblitz<sup>®</sup> CS35

35mm Uni-Stable Optical Shutter

#### Overview

The Uniblitz CS35 has been designed to provide accurate, repeatable exposures for a wide variety of applications including microscopy, video imaging, and more. The slim form-factor provides a 35mm aperture that can be inserted into a 3.33 inch diameter housing. The CS35 is available in a housed or an un-housed configuration for OEM applications (or simply where spatial limitations are a consideration).

Uni-stable shutter devices, like the CS35, require power to hold the blades in the open state.

#### Need Support? Please visit our website or email us at info@uniblitz.com.

Tel: <u>585-385-5930</u> | Toll-Free: <u>800-828-6972</u> | Fax: 585-385-6004 | 803 Linden Ave. Rochester, NY 14625 Updated 2/18 | Datasheet Version 5.2 | ©2018 Vincent Associates

## **Key Features**

- 35mm aperture
- Uni-stable operation
- **<u>501A-S7</u>** adapter included
- Slim, small form-factor
- RoHS Compliant
- Transfer time on opening:

#### 13.0 milliseconds

• Total opening time:

#### 17.0 milliseconds


Configured for the <u>VCM-D1</u>
 <u>Shutter Driver</u>

## **Product Options**

CS35 2 3 4 5 6 -	7 - 8 Ex: CS35S3T0-EC-21
<ul><li>Shutter Series:</li><li>CS35</li></ul>	<ul> <li>Driver Compatibility:</li> <li>S: Use with VCM-D1 (Std.)</li> <li>E: Use with D880C or VED24</li> </ul>
<ul> <li>3 Housing:</li> <li>1: Un-Housed</li> <li>3: #3 Housing</li> </ul>	<ul> <li>4 Blade Coating: <sup>1</sup></li> <li>T: Low Energy (Teflon<sup>®</sup>)</li> <li>ZM: High Energy (AlMgF2) <sup>2</sup></li> </ul>
<ul> <li>Electronic Sync:</li> <li>0: Omitted</li> <li>1: Included</li> </ul>	<ul> <li>Connector:</li> <li>L: 18" flying leads (Un-housed only)</li> <li>Leave blank for 5-pin Switchcraft connector</li> </ul>
<ul> <li>Encapsulated Coil:</li> <li>EC: Included</li> <li>Leave blank if not required</li> </ul>	<ul> <li>Mount: (#3 housing only)</li> <li>21: Zeiss Axiovert</li> <li>24: Olympus BX/IX</li> <li>27: Nikon</li> <li>105: C-Mount (Male)</li> <li>106: C-Mount (Female)</li> <li>Leave blank if not required</li> </ul>

<sup>1</sup> Other blade coating options may be available by special order.
<sup>2</sup> Input side only; Teflon<sup>®</sup> coating is on opposite side to protect shutter blade surface. Light source must be input to the reflective side only.

## Shutter Timing

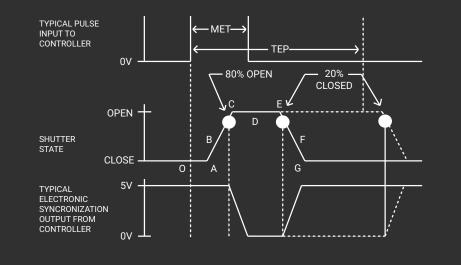

# CS35 (w/ VCM-D1 driver and Teflon® coated blades)Time (msec.)O - ADelay time on opening after current applied4.0A - CTransfer time on opening13.0O - CTotal opening time17.0

A - C	Transfer time on opening	13.0
O - C	Total opening time	17.0
C - E	Min. dwell time with min. input pulse	10.0
B - F	Min. equivalent exp. time	17.5
E - G	Transfer time on closing	30.0
A - G	Total window time	53.0
MET	Min. exposure time	20.0
TEP	Typical exposure pulse	>20.0


#### **Technical Specifications**

Coil	Voltage to	Hold Voltage	
Resistance	Open	(Nominal) <sup>1</sup>	
12 Ω	+70 VDC	+7 VDC / +5 VDC <sup>2</sup>	

<sup>1</sup> Voltage level required across actuator coil when held in the open position.
<sup>2</sup> Dual hold voltage system included in <u>VCM-D1 Driver</u>


<sup>3</sup> (Continuous/Burst) Continuous frequency rating specified at shutter's minimum exposure pulse. Burst frequency rating specified for four (4) seconds maximum with one (1) minute minimum between bursts.

Series	Weight	Weight	Operating	Max. Opening	Max. Closing	Max. Freq. of	Number of
	(Unhoused)	(Housed)	Temp.	Bounce	Bounce	Operation <sup>3</sup>	Shutter Blades
CS35	1.64 oz (0.05 kg)	6.70 oz (0.19 kg)	0 - 80 °C	15%	5%	2.5 Hz / 15 Hz	5

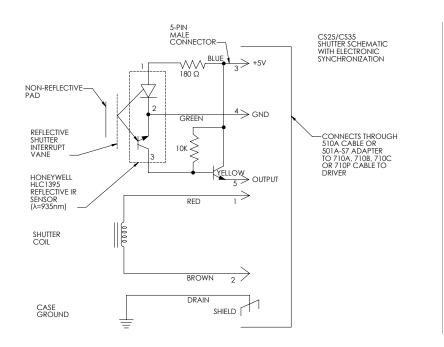


The electronic synchronization system provides a feedback signal (through the driver utilized) after the shutter transfers to the open state. The system incorporates an infrared emitting diode, an infrared sensitive detecting transistor, and an interrupting vane. The vane is attached to the shutter so as to block the light path between the emitter and detector in the closed position. When the shutter transfers to the 80% open position, the vane is removed from the infrared light path, allowing the emitter to switch the detector to the active state. For the CS35, this system uses a similar activation flag attached to the mechanism, which triggers a reflective emitter/detector device. **No connection to the designated synchronization pins when no electronic sync. is selected.** 

#### Technical Drawings - Un-housed CS35

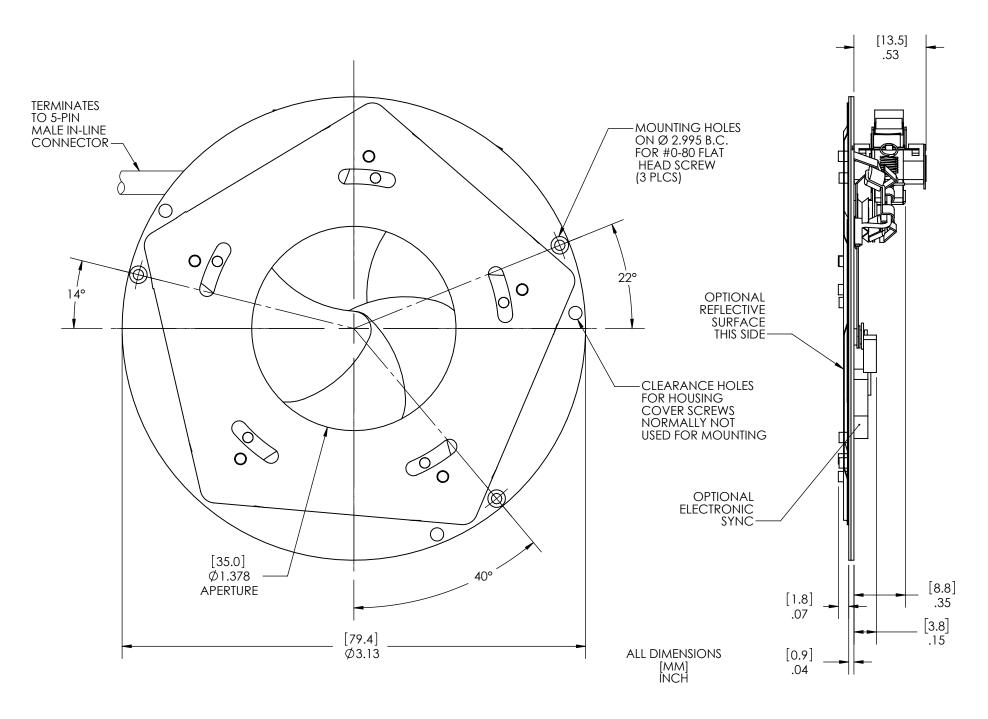

## Technical Drawings - Housed CS35

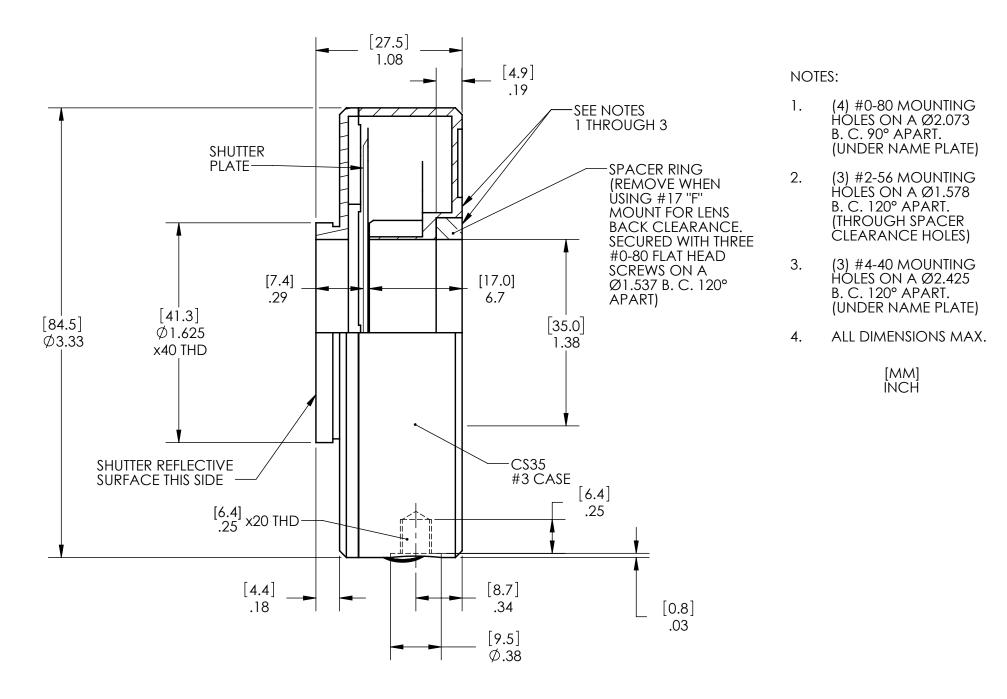

#### Technical Drawings - CS35 Connector Layout

